

## **ATL-SP, Side Ported Compatibility Vessel Instructions**

**THIS EQUIPMENT IS ONLY TO BE USED BY FULLY CERTIFIED  
AND TRAINED PERSONAL. THE USER IS FULLY RESPONSIBLE  
FOR ANY INJURY OR DAMAGE RESULTING  
FROM THE USE OF THIS EQUIPMENT**

Read & follow the safety sheet instructions. In order to fill & use you must remove the vessel from the shield. First re- move the 3 thumb screws from the bottom of the shield. Grab the aluminum bottom part through the slot and re- move it. Now you can remove the vessel & have access to the coupling. When loosening or tightening the coupling ALWAYS remove it from the shield. When installing any hardware to the side port NEVER apply excess force to the threaded stem. Be very careful when washing the vessel. Do not scratch the PVC coating. If the coating is scratched or removed in anyway send it back to us to be recoated. If cracks in the shield appear, have it replaced. If there are any questions or concerns please contact:

Aero-Tech Laboratory Equipment Company, LLC

Phone: (870) 420-3613

Email: [info@aerotechlab.com](mailto:info@aerotechlab.com) Web site: [www.aerotechlab.com](http://www.aerotechlab.com)

Disclaimer:

Aero-Tech Laboratory Equipment Company, LLC

shall not be liable for any injury, loss or damage, direct or consequential, arising out of the use of this product. Due to the nature or research and development, the user is responsible for determining the suitability of the product for it's intended use.