

# Instructions For Aerosol Can Puncturing Device

Make sure the puncturing needle is back up into the valve, so it does not interfere when attaching it to the can. Attach the AC-PD to the can by slipping the inner bottom lip just under the outside curl edge of the valve. Hold it against the valve while holding the can with your thumb or index finger which ever is more comfortable for you. Screw the handle in clockwise until you feel contact, go 2 or more turns. The can valve is now pierced. When you back it out you will be relieving the pressure in the can. If you desire you can screw the handle back in and the leakage will stop. If you do not hold the can straight up the product will leak out. Depending on the type of propellant being used it may be necessary to use an explosion proof hood. All safety precautions must be exercised due to the nature of propellants.

If you have any questions or need parts or service, please call (417) 533-5552 or email us [info@aerotechlab.com](mailto:info@aerotechlab.com)

Aero-Tech Laboratory Equipment Company, LLC  
21397 Tampa Drive, Lebanon, MO 65536 U.S.A.

**Disclaimer:** Aero-Tech Laboratory Equipment Company, LLC shall not be liable for any injury, loss or damage, direct or consequential, arising out of the use of this product. Due to the nature of research and development the user is responsible for determining the suitability of the product for its intended use.