

Instructions For Aerosol Can Puncturing Device

THIS EQUIPMENT IS ONLY TO BE USED BY FULLY CERTIFIED AND TRAINED PERSONAL. THE USER IS FULLY RESPONSIBLE FOR ANY INJURY OR DAMAGE RESULTING FROM THE USE OF THIS EQUIPMENT

Make sure the puncturing needle is back up into the valve, so it does not interfere when attaching it to the can. Attach the AC-PD to the can by slipping the inner bottom lip just under the outside curl edge of the valve. Hold it against the valve while holding the can with your thumb or index finger which ever is more comfortable for you. Screw the handle in clockwise until you feel contact, go 2 or more turns. The can valve is now pierced. When you back it out you will be relieving the pressure in the can. If you desire you can screw the handle back in and the leakage will stop. If you do not hold the can straight up the product will leak out. Depending on the type of propellant being used it may be necessary to use a explosion proof hood. All safety precautions must be exercised due to the nature of propellants.

If you have any questions or need parts or service, please call (870) 420-3613 or email us info@aerotechlab.com

Disclaimer: Aero-Tech Laboratory Equipment Company, LLC shall not be liable for any injury, loss or damage, direct or consequential, arising out of the use of this product. Due to the nature of research and development the user is responsible for determining the suitability of the product for it's intended use.